

CARSON®

MicroMini™

20x Pocket Microscope



MM-280B MM-280G MM-280O

Congratulations on selecting your new MicroMini™ pocket microscope! In order to achieve optimum performance, please follow instructions for proper use and care.

INSTALLING THE BATTERY:

Loosen battery door screw with a screwdriver and remove. (Fig. 1).

Install 3 LR43 batteries with the correct polarity (Fig. 2). Replace battery door and tighten screw with screwdriver.

WARNING:

Use only alkaline batteries. Mixing old and new, different brands or types within brands may cause batteries to leak, overheat or explode. Keep away from fire or excessive heat. Inspect batteries for corrosion or leakage. Insert in direction indicated. Rechargeable batteries not recommended.

USING THE MICROSCOPE:

Place the microscope directly on top of the viewing object and rotate the focusing ring until the image is sharp and clear (Fig. 3).

Slide LED switch (Fig. 4):

- Down to activate the microscope light
- Up to use the microscope as a flashlight
- Turn off LED by sliding to center position

USING THE UV LIGHT:




Press and hold UV light switch to activate UV light (Fig. 5).

CARE AND MAINTENANCE:

Clean using microfiber lens cloth. Do not use any abrasives, chemical cleaners, or solvents.

CAUTION to avoid possible permanent eye damage:

- Do not look at the sun or bright objects with this microscope
- Do not place the lighted microscope directly upon human eye

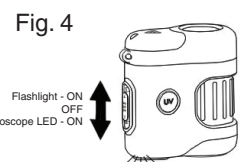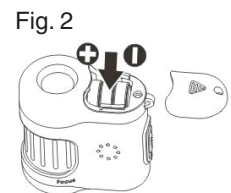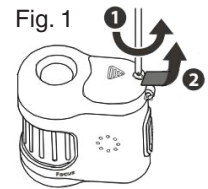

Customer Service:

If you experience any difficulties, please contact us and we will be happy to help you.

In the U.S.A.

Phone: 1-631-963-5000
Email: info@carson.com

In the U.K.

Phone: +44 (0) 1938 557440
Email: uksupport@carson.com

In the E.U.

Phone: +44 (0) 1938 557440
Email: eusupport@carson.com

For warranty information, visit www.carson.com/warranty