
**MIKROSKOP
MICROSCOPE
40X-1280X**



- DE** Schnellstart-Anleitung
- EN** Quick-start instructions
- FR** Guide de démarrage rapide
- NL** Snelstart-handleiding
- IT** Guida rapida
- ES** Guía de inicio rápido
- RU** Краткое руководство пользователя





Lieferumfang · Contents · Contenu de la livraison · Inhoud · Accessori in dotazione
Artículos incluidos en la entrega · Material fornecido · Комплект поставки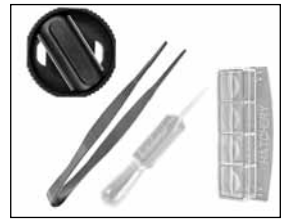

Zubehör anbringen · Accessory installation · Fixation des accessoires · Accessoires aansluiten · Montaggio degli accessori · Instalación de accesorios · Установка дополнительного оборудования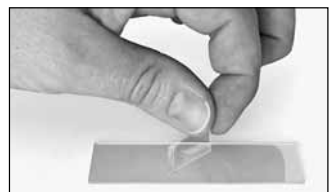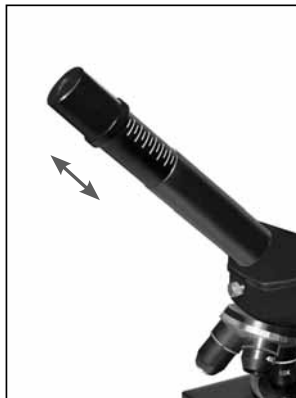

Stromversorgung / Beleuchtung · Power supply and illumination operation · Alimentation électrique / Eclairage · Stroomvoorziening / Belichting · Alimentazione elettrica / Illuminazione Alimentación de corriente / Iluminación · Питание от сети / Подсветка

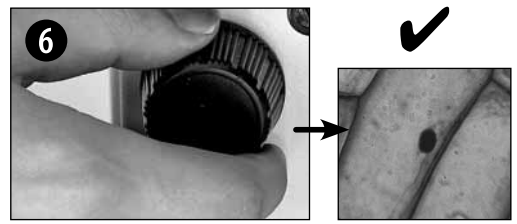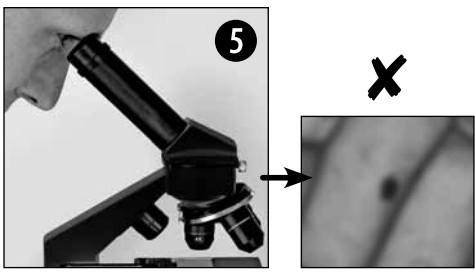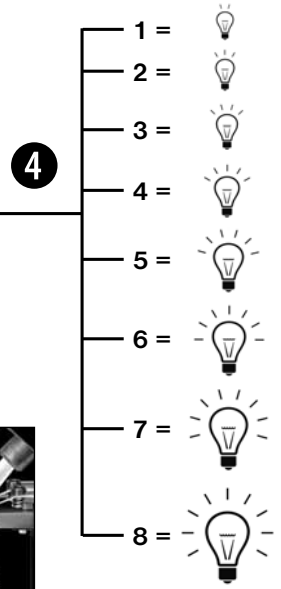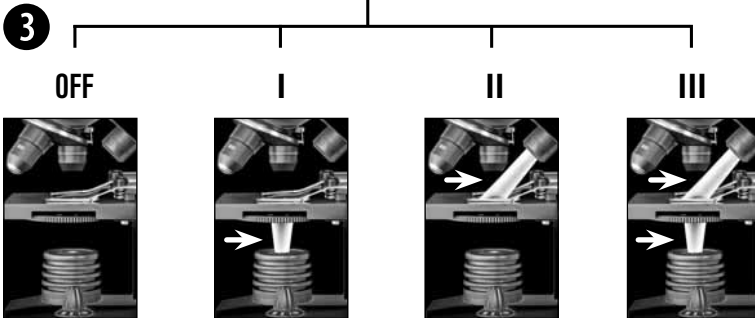Download: <http://www.bresser.de/9039000>

Eine ausführliche Anleitung steht für Sie auf unserer Internetseite zu Ihrer Verfügung.
A more detailed manual is available on our website. • Un guide détaillé est disponible sur notre site. • Een uitgebreide handleiding is beschikbaar op onze website. • Un manuale di istruzioni dettagliato è disponibile sul nostro sito Web su internet. • Un manual de Instrucciones detallado se encuentra disponible en nuestra página Web en internet.
Подробное руководство можно найти на веб-сайте.

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EN About this manual

This quick-start guide will help you put the device into operation quickly, but it is not intended to replace the more detailed user manual.

Please read the safety instructions and the user manual carefully before using the device. This quick-start manual contains the most important steps to start using your new device. Comprehensive operating instructions and further information, including troubleshooting advice, is available on our website: www.bresser.de (Download: <http://www.bresser.de/9039000>)

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General Warning

DANGER!

Tools with sharp edges and points are often used when working with this device. Because there is a risk of injury from such tools, store this device and all tools and accessories in a location that is out of the reach of children.

This device contains electronic components which operate via a power source (power supply). Do not let children use the device while unattended. Use the device only as described in the manual, otherwise you run the risk of an electric shock.

Do not expose the device to high temperatures. Use only the power adapter supplied. Do not short-circuit the device or batteries, or throw them into a fire. Excessive heat or improper handling could trigger a short-circuit, a fire or an explosion.

Never bend, pinch, pull or run over the power and connecting cables, extensions and adapters. Protect the cables from sharp edges and heat.

Before operating, check the device, cables and connections for damage. Never use a damaged unit or a unit with damaged power cables. Damaged parts must be exchanged immediately by an authorised service centre.

Children should only use the device under adult supervision. Keep packaging material, like plastic bags and rubber bands, out of the reach of children, as they pose a choking hazard.

CAUTION!

Children must not have access to the included chemicals and liquids. Do not drink the chemicals. Wash hands thoroughly with running water after using the chemicals. In the event that the chemicals come into contact with your eyes or mouth, rinse thoroughly with water. If you are in pain after exposure, contact a doctor immediately and take the substances with which you came into contact with you.

Before cleaning the device, disconnect it from the power supply by removing the plug.

NOTE!

Position your device so that it can be disconnected from the power supply at any time. The wall socket you use should be located near the device and easily accessible, since the plug on the power cable serves as a disconnecting device for the power supply.

Always pull on the plug to separate the device from the power supply. Never pull on the power cable.

Do not disassemble the device. In the event of a defect, please contact your dealer. The dealer will contact the Service Centre and can send the device in to be repaired, if necessary.

Do not expose the device to temperatures above 45 °C.

Declaration of Conformity

CE Bresser GmbH has issued a 'Declaration of Conformity' in accordance with applicable guidelines and corresponding standards. This can be viewed any time upon request.

Warranty and warranty term extension

The warranty term is two years from the date of purchase. Please keep your proof of purchase. Register at **www.bresser.de/warranty** and fill out a brief questionnaire to get your warranty term extended to **five years**. Registration must be completed within three months of purchase (date of receipt) to validate the warranty. If you register thereafter, the warranty term will not be extended.

If you have problems with your device, please contact our customer service first. Do not send any products without consulting us first by telephone. Many problems with your device can be solved over the phone. If the problem cannot be resolved by phone, we will take care of transporting your device to be repaired. If the problem occurred after the warranty ended or it is not covered by our warranty terms, you will receive an estimate of repair costs, free of charge.

Service Hotline: +49 (0) 2872 - 80 74-210

Important for any returns:

Please make sure to return the device carefully packed in the original packaging in order to prevent damage during transport. Also, please enclose your receipt for the device (or a copy) and a description of the defect. This warranty does not imply any restriction of your statutory rights.

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are trademarks of the National Geographic Society,
used under license. All rights reserved.